

Milk Security Test

Detection of
NEUTRALIZERS
HYDROGEN PEROXIDE
UREA

in raw milk

12 tubes / 600
strip tests


**Rapid & sensitive
dry strips to detect
milk adulteration**

**Pure milk
for
the Dairy industry**


Test in three simple steps:

STEP 1

Place the test strip into the milk sample.
Attention! The milk must cover the indication areas.

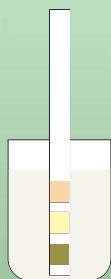

STEP 2

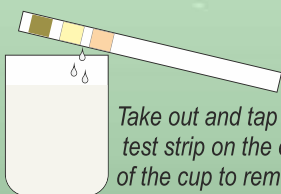

Take out and tap the test strip on the edge of the cup to remove excess milk.

Wait
60 seconds



STEP 3

Compare the indication areas of the test strip with the color scale on the label.

NEUTRALIZERS (%) 120s	0	0.03	0.06	≥0,1
	0	0.0075	0.0015	≥0,025
HYDROGEN PEROXIDE (%) 80s	0	0.07	≥0,12	
	0	0.07	≥0,12	
UREA & NEUTRALIZERS (%) 90s	0	0.07	≥0,12	
	0	0.07	≥0,12	

**Box dimensions
(W x D x H):
115 x 155 x 100 mm
Weight < 0.5 kg**

Carton storage box



12 tubes / 600 strip tests

**Rapid & sensitive
dry strips to
determine
titratable acidity**

**High-quality milk
for
the Dairy industry**

**Box dimensions
(W x D x H):
115 x 155 x 100 mm**

Weight < 0.5 kg

Carton storage box

**buy online on
www.lactoscan.com**



Milk Security Test

Determination of
**TITRATABLE
ACIDITY**

in raw milk




Test in three simple steps:

STEP 1

Place the test strip into the milk sample.
Attention! The milk must cover the indication areas.

STEP 2

Take out and tap the test strip on the edge
of the cup to remove excess milk.

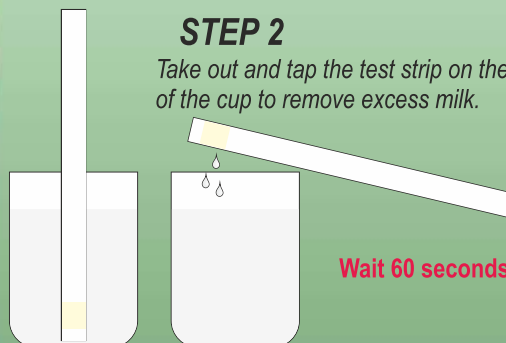

Wait 60 seconds

STEP 3

Compare the indication areas of the
test strip with the color scale on the label.

15,0	18,0	24,0	27,5	Thörner degrees, °Th
13,5	16,2	21,6	24,75	Dormic degrees, °D
6,0	7,2	9,6	11,00	Soxhlet Henkel degrees, °SH
6,9	6,6	6,3	6,00	pH

MILKOTRONIC LTD

Bulgaria, 6000 Stara Zagora
167, Tsar Simeon Veliki, A, 25
Tel: + 359 457 670 82,
Fax: + 359 457 670 83

e-mail: salesdpt@lactoscan.com
Eng, Spa, Port: +359 886 841 389
e-mail: sales2@lactoscan.com
Eng, Rus: +359 886 841 395
e-mail: mainservice@lactoscan.com
Eng, Rus: +359 888 811 017

e-mail: m.mihneva@lactoscan.com
Eng: + 359 885 597 232
e-mail: milkanalyzers@lactoscan.com
Eng, Rus: +359 896 753 803

distributors: